



Dear customer

You have received an offer from us for a machine with natural refrigerant (NR) or placed an order for such a machine.

In the standards and regulations are rules concerning machines containing natural refrigerants. In this letter, we wish to make you aware of these particular conditions. In addition we wish to inform you concerning attention to operation within these regulations and to operate within the guidelines set out in the relevant operating manual.

- 1) The machine was built in accordance to the EU and the EFTA states regulations.
- 2) Setting up the machine in public buildings is not allowed.
- 3) The machine is not certified for use in an ATEX area.
- 4) Also for smaller machines it is necessary that the size of the operating environment and the quantity of refrigerant must be matched. The size of the room required is calculated as follows:

**max. 8g Refrigerant / m<sup>3</sup> Room size.**

Example:

A machine with 100 g. refrigerant has to be placed in a room bigger then 5 m<sup>2</sup> floor space if the height is 2,5m.

In reality, this will practically never happen, but still we want to support you to keep your compliance with the effective standards.

Take the quantity of refrigerant used from the data sheet or the type number label on the rear of the machine. Data sheets can be obtained via the internet at

[www.huber-online.com](http://www.huber-online.com)