Inspired by **temperature**

Software FlowHow+

Handbuch

huber

Software FlowHow+

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V1.0.0de/19.05.21//V2.1.0

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2.1 Thermofluid manuell auswählen

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The interface consists of a top section with 'Load *.fcv' and 'Save *.fcv' buttons. Below is a table with two columns: 'Temperature in °C' and 'Viscosity in mm²/s'. The table has 14 rows. Below the table are radio buttons for 'Viscosity' and 'Density', and buttons for 'Graph', 'Clear', and 'Activate'. A 'Fluid:' text box is also present.

Left Screenshot (Initial State):

- Buttons: Load *.fcv, Save *.fcv
- Table headers: Temperature in °C, Viscosity in mm²/s
- Table rows: 1 to 14 (empty)
- Buttons: Viscosity, Density, Graph, Clear
- Fluid: (empty text box)
- Buttons: Activate
- Annotations: [A] points to Temperature header, [B] points to Viscosity header, [C] points to Save *.fcv, [D] points to Fluid text box, [E] points to Activate button.

Right Screenshot (Activated State):

- Buttons: Load *.fcv, Save *.fcv
- Table headers: Temperature in °C, Viscosity in mm²/s
- Table data:


	Temperature in °C	Viscosity in mm²/s
1	-50	6
2	-40	5.2
3	-30	4.3
4	-20	3.5
5	-10	2.8
6	0	2.2
7	10	1.8
8	20	1.5
9	30	1.3
10	40	1.1
11	50	0.95
12		
13		
14		
15		
- Buttons: Viscosity, Density, Graph, Clear
- Fluid: Ethanol
- Annotation: [F] points to the right side of the table.

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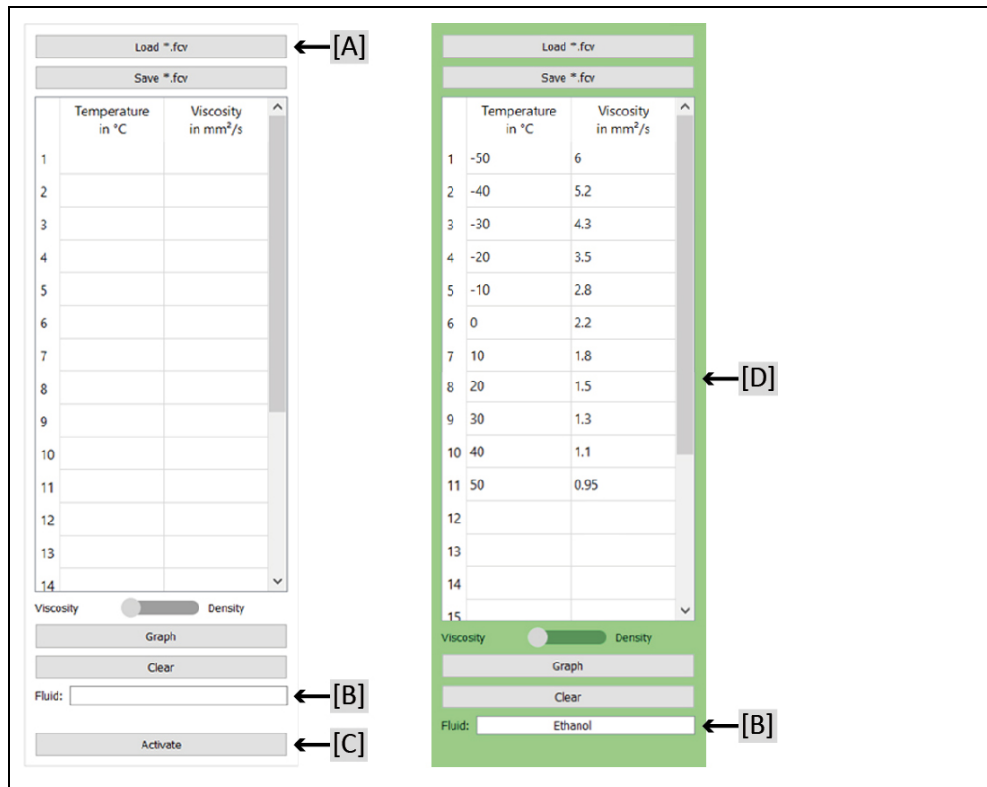The interface consists of two main panels. The left panel is the initial state, and the right panel shows the state after loading a file.

Left Panel (Initial State):

- Buttons: Load *.fcv [A], Save *.fcv
- Table with 14 rows and 2 columns: Temperature in °C, Viscosity in mm²/s
- Controls: Viscosity/Density toggle, Graph, Clear, Fluid: [B], Activate [C]

Right Panel (Loaded State):

- Buttons: Load *.fcv, Save *.fcv
- Table with 15 rows and 2 columns: Temperature in °C, Viscosity in mm²/s
- Controls: Viscosity/Density toggle, Graph, Clear, Fluid: Ethanol [B]
- The entire right panel is highlighted with a green border, and the data table is labeled [D].

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Inspired by **temperature** designed for you

Peter Huber Kältemaschinenbau AG
Werner-von-Siemens-Str. 1
77656 Offenburg / Germany

Telefon +49 (0)781 9603-0
Telefax +49 (0)781 57211

info@huber-online.com
www.huber-online.com

Technischer Service: +49 (0)781 9603-244

-125 °C ... +425 °C

huber